

ELVATECH PRECISION THAT
EMPOWERS CONFIDENCE

ELVAX PLUS



**MEET THE NEW XRF STANDARD WITH
THE NEWEST BENCHTOP X-RAY
FLUORESCENCE SPECTROMETER**

ELVAX PLUS SPECTROMETER IS A DESKTOP ENERGY-DISPERSIVE X-RAY FLUORESCENCE (EDXRF) ANALYZER WITH EXTENDED ELEMENT RANGE FROM NA (Z=11) TO U (Z=92).

ElvaX Plus spectrometer is a desktop energy-dispersive X-ray fluorescence (EDXRF) analyzer with extended element range from Na (Z=11) to U (Z=92). The Helium purge system and the automatic primary filter changer combined with the large area SDD detector are used to allow effective registration of the light elements spectra.

- ✓ World's fastest XRF analyzer (throughput above 500 000 cps) – highest accuracy and excellent precision within the shortest measurement time;
- ✓ Wide element detection range from Na (Z=11) to U (Z=92);
- ✓ Helium purge feature improves sensitivity to light elements (Na, Mg, Al, Si, P, S);
- ✓ Excellent calibration stability, compensation of ambient temperature and pressure variations;
- ✓ Optional 16-position automatic sample changer increases productivity of your lab;
- ✓ Optional high-resolution CCD camera for precise sample positioning;
- ✓ Low maintenance cost.



ACCURACY

With innovative hardware and unique software algorithms, the ElvaX Plus offers the highest precision and accuracy of elemental analysis for a whole range of analytical applications.

SPEED

Dynamically Adaptive Shaping (DAS) DPP developed by Elvatech combined with the ultra large area Fast SDD detector makes the ElvaX Plus the fastest XRF analyzer on the market.

STABILITY

Thanks to the **digiX-50** digital X-Ray source, automatic temperature and pressure corrections, and automatic calibration adjustment, ElvaX Plus offers long-life reproducibility of the analysis without the need to re-calibrate the system.

ElvaX Plus spectrometer is used for quantitative and qualitative analysis of the elemental composition of metal alloys, liquids, powders, food and biological samples, as well as samples deposited on surfaces or filters, in a wide range of element concentrations.

Analyzer's detecting limit is better than 1 ppm for most elements in a light matrix.



The optional automatic 16-positions sample changer coupled with the fast measurement increases the productivity of your lab up to 1000 samples per workday.

At. Numb	Element	Series	Intensity	Concentration
28	Ni	K	145153	63.2316 ± 0.2404%
27	Co	K	18370	8.6076 ± 0.1553%
42	Mo	K	40521	6.4280 ± 0.0692%
74	W	L	6766	6.3900 ± 0.4055%
24	Cr	K	8804	6.2659 ± 0.1517%
26	Fe	K	10730	4.3035 ± 0.0562%
22	Ti	K	611	1.3837 ± 0.2246%
52	Te	L	3688	0.8797 ± 0.0804%
17	Cl	K	1891	0.8478 ± 0.2206%
19	K	K	2975	0.4896 ± 0.0610%
23	V	K	412	0.4766 ± 0.1416%
79	Au	L	594	0.3176 ± 0.1564%
25	Mn	K	574	0.2301 ± 0.1463%
13	Al	K	824	0.0822 ± 0.0076%
14	Si	K	868	0.0232 ± 0.0070%

The superior throughput, which is provided by state of the art DAS DPP, makes it possible to get the highest accuracy analysis results in just a few seconds.

DIGITAL X-RAY SOURCE DIGIX-50

Anode: W, Rh, Ag

Voltage: 50 kV

Current: 200 μ Amp

Power: 5 W

5 position filter changer

X-RAY DETECTOR

Type: Large Area SDD

Area: 20 mm²

Energy resolution: < 140 eV at Mn K α ,
< 85 eV at Al K α

Count rate: 500 000 cps

ELECTRONICS

DPP: proprietary DAS (Dynamically Adaptive Shaping) type, 80 MHz sampling rate

MCA: 4096 channels

GENERAL

Dimensions: 430 x 340 x 200 mm

Weight: 18 kg

Power supply: 90 – 240 V, 50/60 Hz

Power consumption: 50 W

SOFTWARE

Operating system: Windows XP/Vista/7/8/10

Analysis algorithms: Fundamental parameters (FPA), Empirical (regression) calibrations, Manual spectra comparison



Elvatech Ltd. 50, Mashynobudivna str., Kyiv, 03067, Ukraine

Phone: (+1929) 243 99 06, (+38 063) 576 21 21, (+38 095) 576 21 21,

E-mail: office@elvatech.com

www.elvatech.com

